**Modular Series**

# INSTRUCTION MANUAL

Model
SZ-STL1
SZ-STL1H
SZ-STL2
SZ-STL2H

v 1.0 2018

---

## **SUMMARY**

**Warning**

**Intended use**

**Overview**

**Assembly**

**Using the base**

**Equipment disposal**

## Warning

This microscope is a scientific precision instrument designed to last for many years with a minimum of maintenance. It is built to high optical and mechanical standards and to withstand daily use.


We remind you that this manual contains important information on safety and maintenance, and that it must therefore be made accessible to the instrument users.

We decline any responsibility deriving from incorrect instrument use uses that does not comply with this manual.


## Intended use

For research and teaching use only. Not intended for any animal or human therapeutic or diagnostic use.

## Overview SZ-STL1 and SZ-STL1H

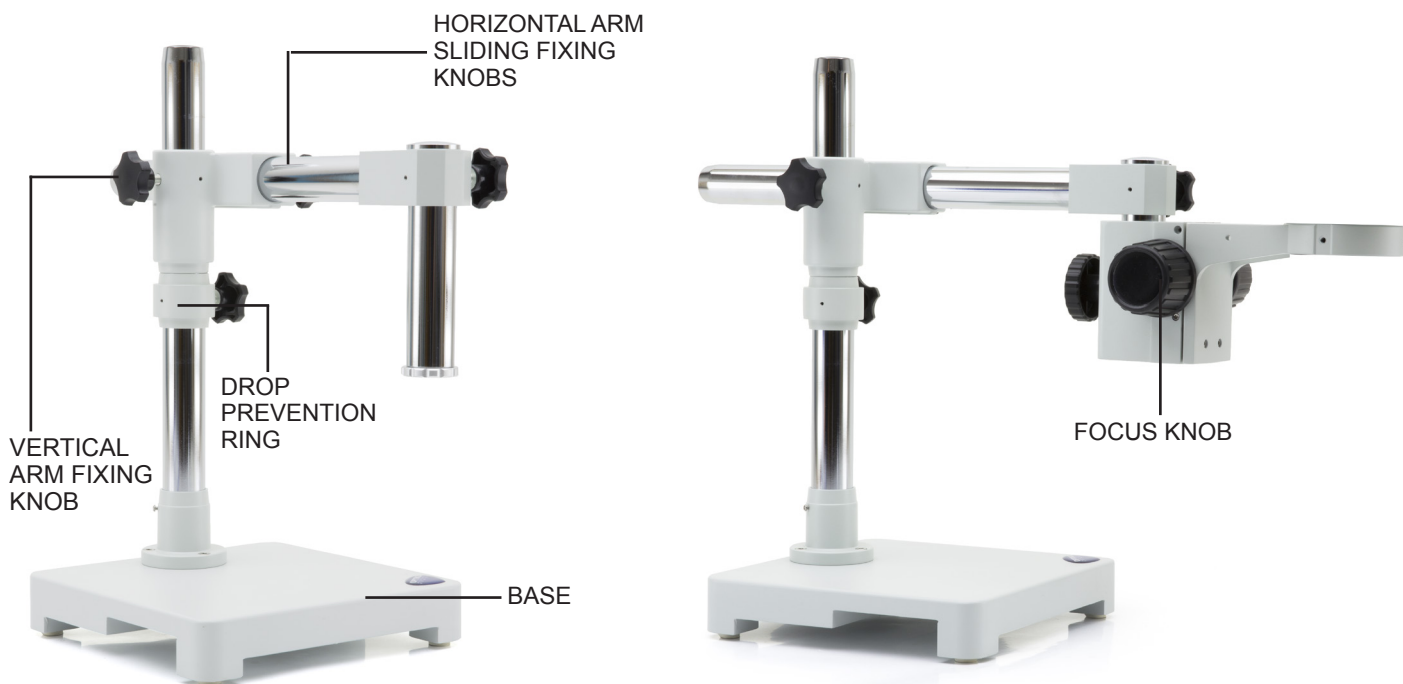## Overview SZ-STL2 and SZ-STL2H
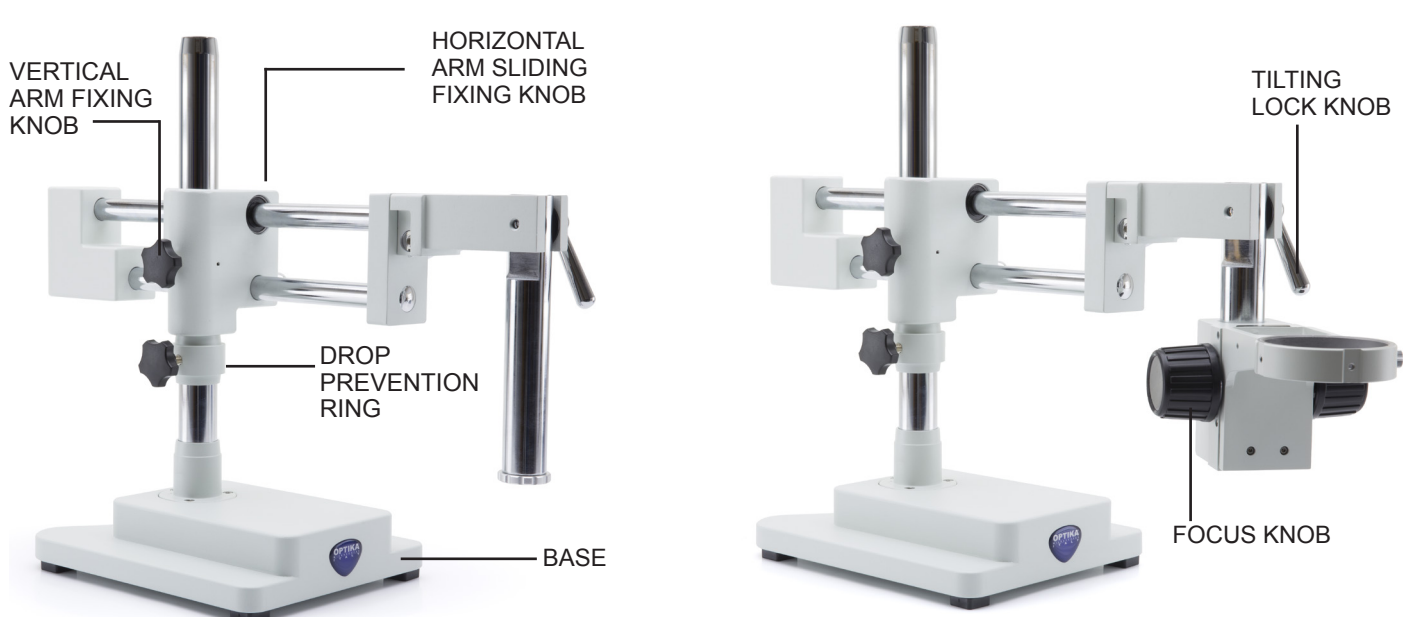
## Assembly (SZ-STL1 / SZ-STL1H)

1. Screw the pillar on the base. (Fig. 1)
2. Tighten the screw to lock the pillar. (Fig. 2)
3. Insert the drop preventing ring. (Fig. 3)
4. Insert the horizontal arm and lock it with the locking knob. (Fig. 4 & 5)
5. Install the focus holder. Unscrew the prevention knob ②, and insert the head holder arm ① from above in the hole of the horizontal arm. (Fig. 6-7)

► **ONLY FOR SZ-STL1H:**

6. Install the head holder, lock the fixing knob and screw the prevention knob ② from below. (fig. 8-9)

## Assembly (SZ-STL2 / SZ-STL2H)

1. Screw the pillar on the base (Fig. 10)
2. Tighten the screw to lock the pillar (Fig. 11)
3. Insert the drop preventing ring. (Fig. 12)
4. Insert the horizontal arm and lock it with the locking knob ① (Fig.13). For a bigger safety, lock the fixing screw ② with the provided allen wrench (Fig. 14)

► **ONLY FOR SZ-STL2H:**

6. Install the head holder. Unscrew the prevention knob ① and insert the head holder ② from below in the vertical arm. (fig. 15-16). At the end screw again the prevention knob.

## Using the base

### SZ-STL1 / SZ-STL1H:

Moving the horizontal arm (Fig. 17).  
Unlock the knob on the right side of the horizontal arm ①.

The arm can be extended or shortened according to specific needs. (Fig. 18-19)

Fig.17

Fig.18

Fig.19

Swivel the head (Fig. 20)  
Unlock the fixing knob ① and rotate the head to the desired swivel angle (left or right), then fix the knob again.

Fig.20

### Rotating the horizontal arm (Fig. 21)

Loosen the horizontal arm fixing knob ② and rotate the arm, then fix again the fixing knob.

- ▶ **NOTE: 180° rotation of the microscope with respect to the base could cause a rollover of the entire system.**

Fig.21

### Adjusting the tension of the focus knob (Fig. 22)

- ▶ **This adjustment allows to increase or decrease the tension of the knob by avoiding an involuntary descent of the microscope body under its own weight. Adjust the tension just above the point where the focus is stable.**

Grab the knobs ③ with both hands and, holding right knob rotate the left one. Tension increases or decreases according the sense of rotation of the left focus knob.


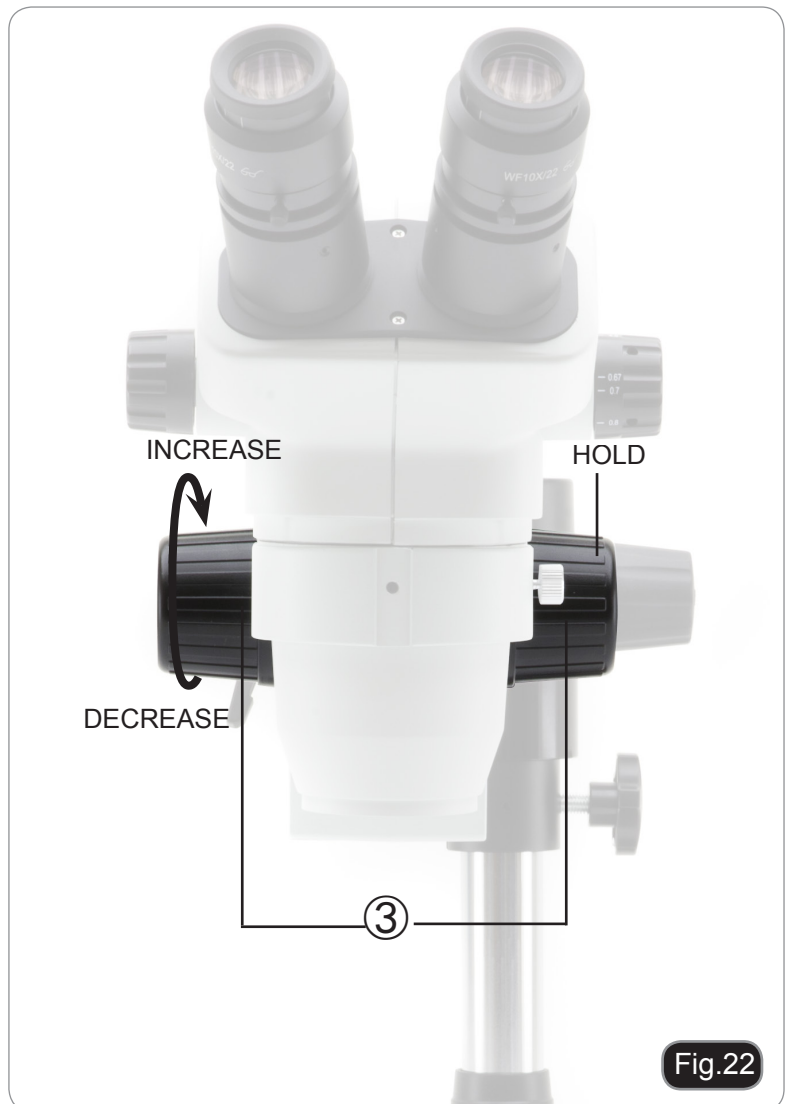
Fig.22

## Using the base

### SZ-STL2 / SZ-STL2H:

Moving the horizontal arm (Fig. 23).  
Unlock the knob on the left side of the horizontal arm ①.

The arm can be extended or shortened according to specific needs. (Fig. 24-25)

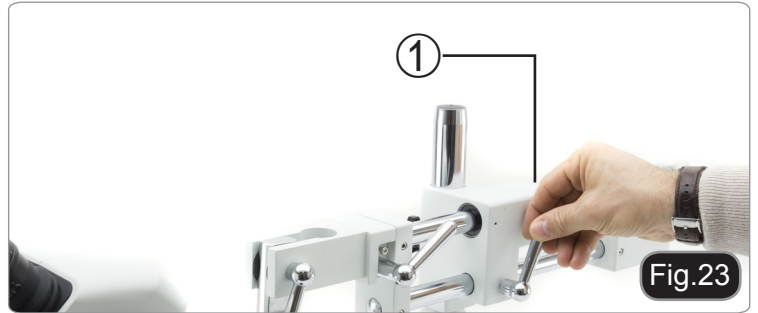Tilting of the head holder. (Fig. 26)  
Loosen the knob ② on the right side of the head holder. The tilting vertical arm can be moved. Once the desired angle is achieved, tighten the wrench again.

Swivel the head (Fig. 27-28)  
Unlock the fixing knob ③ and rotate the head to the desired swivel angle (left or right), then fix the knob again.

Rotating the horizontal arm (Fig. 29)

Loosen the horizontal arm fixing knob ④ and rotate the arm, then fix again the knob.

► **NOTE: 180° rotation of the microscope with respect to the base could cause a rollover of the entire system.**

## Equipment disposal

Art.13 Dlsg 25 July 2005 N°151. "According to directives 2002/95/EC, 2002/96/EC and 2003/108/EC relating to the reduction in the use of hazardous substances in electrical and electronic equipment and waste disposal."

The basket symbol on equipment or on its box indicates that the product at the end of its useful life should be collected separately from other waste.

The separate collection of this equipment at the end of its lifetime is organized and managed by the producer. The user will have to contact the manufacturer and follow the rules that he adopted for end-of-life equipment collection.

The collection of the equipment for recycling, treatment and environmentally compatible disposal, helps to prevent possible adverse effects on the environment and health and promotes reuse and/or recycling of materials of the equipment.

Improper disposal of the product involves the application of administrative penalties as provided by the laws in force.

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**OPTIKA® S.r.l.**